

EMAN Tutorial Session #1

Steve Ludtke

EMAN Documentation

- <> denotes a parameter to fill in, ie <input file>
- [] denotes an optional parameter, ie [mask=<radius>]
- *italics* denote something to be typed into the computer example:

proc2d <input file> <output file> [mask=<r>] [mrc] [spider] [pgm] proc2d file.hed file.spi spider mask=22

- *<program> help* (won't work with python scripts yet)
- <program> (shows usage information)

GUI (eman browser)

- cd data/samples
- run 'eman'

GUI

• v2


• v4

• qplot


- eman
- boxer
- helixboxer
- ctfit
- qsegment(/chimera)
- qindex
- glmatrix

- Non-EMAN Programs
- chimera

The Reconstruction Process

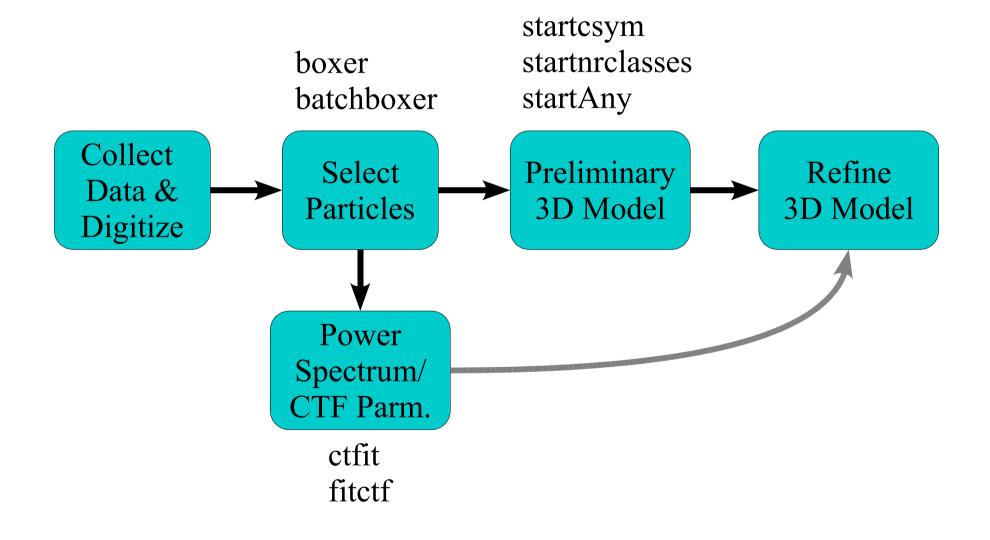

Typical Refinement - EMAN

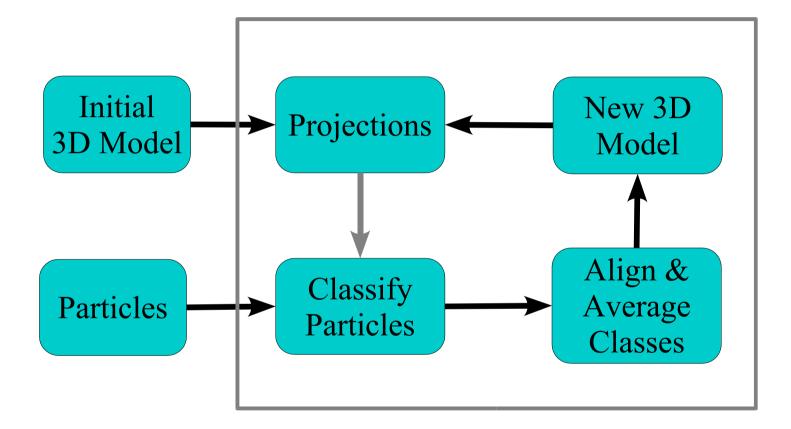

'refine' command

Boxing (aka Particle Selection)

- cd ~/data/GroEL/stage1
- ls
- cat 000script
- boxer (one of the mrc files)

When you get bored ...

• ./000script


CTF Correction


- cd ../stage2
- cat 000script
- ./000script
- ctfit *.tnf
- ./000script2
- eman
- look at files in ../raw and ../raw.w


Initial Model (with symmetry)

- cd ../stage3
- cat 000script
- ./000script
- wait a minute or two
- head start.hed
- head start.filt.hed
- eman
- look at:
- avg.img, ali.img, cls0000.lst, cls0001.lst, classes.img
- sym.img, threed.Oa.mrc, threed.Ob.mrc (final result)

Assume D7 Symmetry

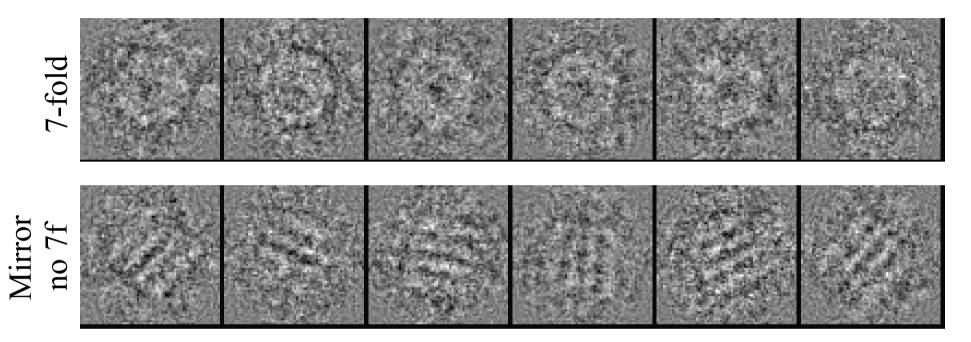

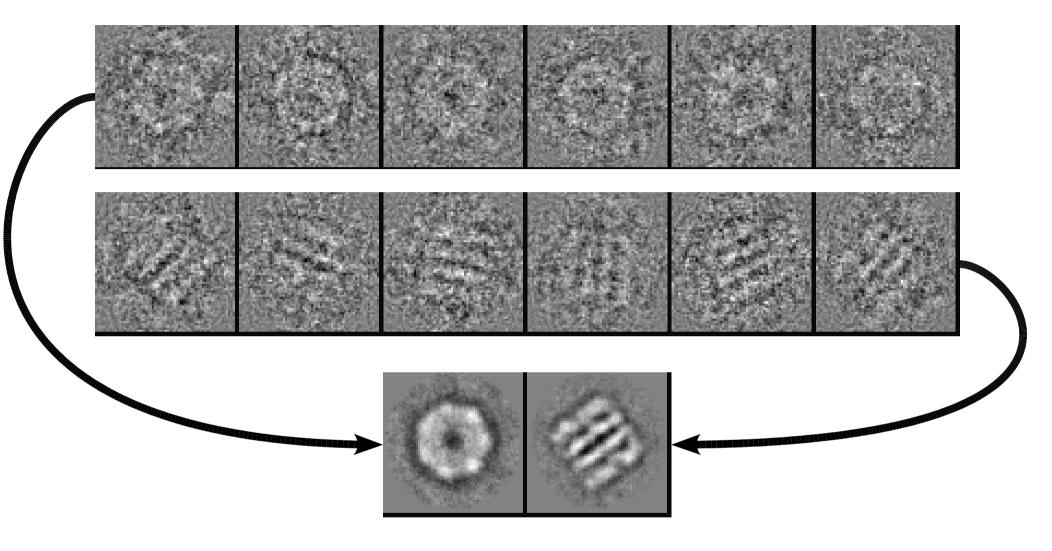

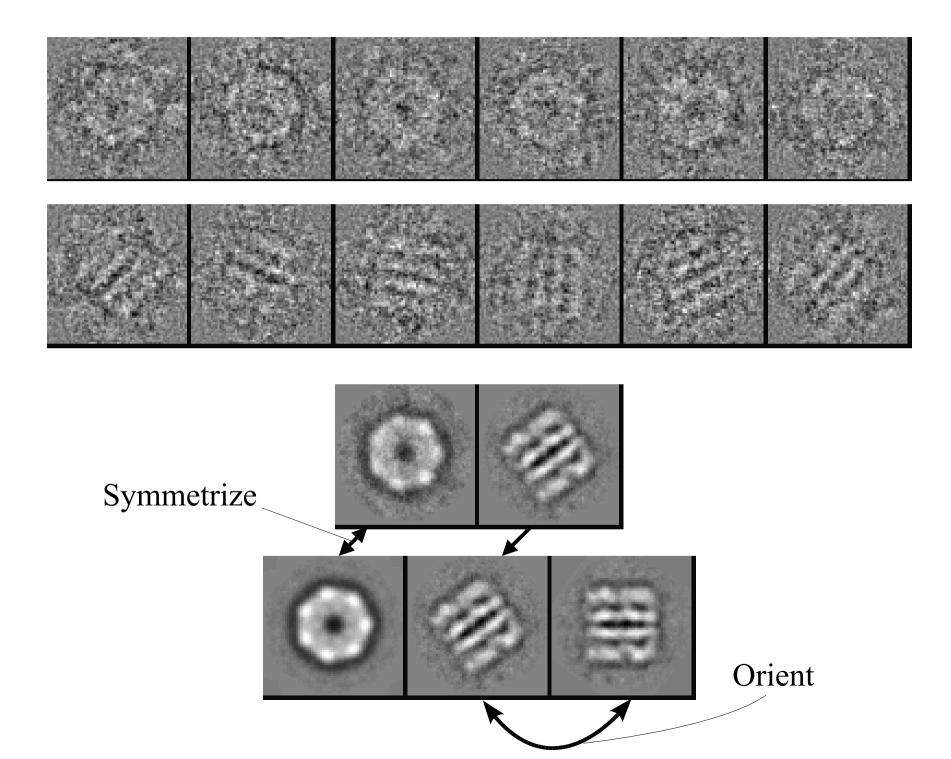

Quick Initial Refinement

- cd ../stage4
- cat 000script
- ./000script
- wait a while

While we wait ...

Boxing (aka Particle Selection)

- cd samples/boxer
- boxer <image>

Tricks:

- *makeboxref.py groel.mrc sym=d7 invert ang=15* (creates a set of reference images from a 3D model for autoboxing)
- batchboxer input=jj5339.f.mrc auto=.3,.7,.1 dbout=5339.box output=5339.ptcl.hed refimg=best.hed

(autoboxes a single micrograph, parameters determined in boxer, refing specifies reference particles for boxing)

Boxing (aka Particle Selection)

- cd /tmp/USERNAME/groel/stage1
- batchboxer input=jj5337.f.mrc dbbox=5337.box output=../stage2/5337.hed invert (takes a list of box locations and chops out each particle into a single image stack)

etc.

(or 'source script', to run the command for all micrographs)

Basic Image Processing

- *iminfo <file> [all] [stat]* (image information)
- proc2d <infile> <outfile> [options] (generic 2d image processing)
- proc3d <infile>
- proc3d <infile> <outfile> [options] (generic 3d image processing)
- procpdb.py <infile> <outfile> [options] (simple PDB file processing)

Quick Initial Refinement

- cd ../stage4
- cat 000script
- ./000script
- wait a while
- eman
- look at:
- classes.*.img, x.*.mrc
- chimera animscript.py

High Resolution Refinement

- cd ../stage5
- cat 000script
- precomputed (looong job)
- eman
- look at:
- classes.*.img, x.*.mrc
- analysis -> convergence
- chimera animscript.py

Simple Docking and Segmentation

- *cd* ../*stage6*
- ./000script
- *chimera fh.0.ent shrunk.mrc ../stage5/threed.8a.mrc*
- ./000script2
- use qsegment
- use chimera

Easier and better ways of doing this will be discussed later in the workshop.

```
• Projection:
mask=<r>
proc=<n>
sym=<cn,dn,etc>
ang=<dang>
tree=<2,3>
perturb
```

Classification: dfilt, phasecls, fscls refine sep=<n> tree=<2,3> shrink=<n> maxshift=<rad> precen, usefilt, slow mask=<r>, imask=<r> amask=<r>, <thr>, <iter> proc=<n>

```
ctfc=<res>, ctfcw=<SF> or median
```

 Class Averages: dfilt, phasecls, fscls refine classiter=<n> classkeep=<sig mult, but not 0> mask=<r>, imask=<r>, imask=<r>, imask=<r>, imask=<r>, othr>, iter> proc=<n> ctfc=<res>, ctfcw=<SF> or median maxshift=<r> precen, slow euler2=<oversmp>

- 3D Reconstruction: hard=<phase err> pad=<size> sym=<cn,dn,etc> 3dit=<n>,3dit2=<n> mask=<r>,imask=<r> proc=<n> collapse=<alt thr>
- Postprocessing (betw iter): xfiles=<A/pix>,<mass in kd>,<ali to> amask=<r>,<thr>,<iter> filt3d=<lp rad>

Typical options to use for a highresolution refinement (EMAN 1.6): mask proc (for clusters) hard ctfcw (requires a SF file) sym (exclude if no symmetry) ang pad classkeep classiter (3 or 0 at the end, 1-2 invalid) xfiles amask (must also use xfiles) refine dfilt (must also use ctfcw) usefilt (start.filt.hed must exist) sep (with oversampled ang=)

REQUIRED:

- <total iter> Final number of iterations you wish to be complete in this directory. ie - if 5 are complete and you specify 6, 1 more iteration will run
- mask=<rad> A circular mask to be applied virtually everywhere. Should be a few pixels larger than the largest radius of your particle
- ctfc=, ctfcw= OR median CTF correction options. ctfc takes a filter resolution in Angstroms, but ctfcw is far superior. ctfcw takes the name of a 1D structure factor file used when fitting the data (MUST be the same file). median does no CTF correction at all. Naturally the data must be properly preprocessed for this option to function, otherwise crashes or invalid results are likely.

- hard=<phase err> Rather obscure. This option determines when a class-average should be excluded from the 3D reconstruction process. 25 is generally good. see 'make3d' for more info.
- sym=<cn,dn,oct,icos> For asymmetric objects, either omit this specification, or use 'c1'. cn denotes a single n-fold rotational symmetry (about the z-axis), dn denotes n-fold dihedral symmetry (cn with n 2-folds in the x-y plane), oct is octahedral (2-3-4, symmetry of a cube), icos is icosahedral (2-3-5).
- ang=<dang> Angular spacing between projections. Smaller numbers produce more projections, usually between 2-10.

REQUIRED:

- pad=<size> This is used to reduce artifacts in Fourier reconstruction. Should be about 25% larger than your model in most cases, and have small prime factors, ie - model size 50 -> pad=64, 70->96, 110->128
- classkeep=<sig mult> This determines how many raw particles are discarded for each class-average. This is defined in terms of the standard-deviation of the self-similarity of the particle set. A value close to 0 (should not be exactly 0) will discard about 50% of the data. 1 is a typical value, and will typically discard ~10-20% of the data.
- refine Almost always specified. This will increase the alignment accuracy (with a speed penalty). May be omitted for early rounds of refinement.

- classiter=<n> Generation of classaverages is an iterative process. Rather than just aligning the raw particles to a reference, they are iteratively aligned to each other to produce a class-average representative of the data, not of the model. ie - this eliminates initial model bias, typically 8 in the early rounds and 3 in later rounds, 0 may be used at the end, but model bias may result.
- dfilt, fscls, phasecls Mutually exclusive options that determine similarity criteria used for classification and quality comparisons. dfilt requires ctfcw, but generally produces the best results. phasecls uses mean phase error. fscls uses Fourier ring correlation. dfilt is an optimized variance with a matched filter.

OPTIONAL:

- filt3d=<rad> Applies a lowpass filter to the 3D model between iterations. This can be used to correct problems that may result in high resolution terms being upweighted. <rad> is the same as for the 'lp=' option in proc3d.
- sep=<n> This interesting option causes each particle to be assigned to the n best classes, not just the single best class. This may be used to smooth and improve fine details in the final stages of a high resolution refinement. Generally used with oversampled ang= values.
- tree=<2,3> This can be a risky option, but it can produce dramatic speedups in the refinement process. Rather than comparing each particle to every reference, this will decimate the reference population to 1/4 or 1/9 of its original size, classify, then locally determine which of the matches is best. Is is safest in conjunction with very small angular steps, ie - large numbers of projections. The safest way to use this is either a) for high-resolution, smallang refinement or b) for the initial iterations of refinement (then turn it off for the last couple of iterations). May not work on certain cluster configurations.

TYPICAL (not required):

- xfiles=<A/pix>,<mass>,<ali to> This is a convenience option. For each 3D model it will produce a corresponding x-file: threed.1a.mrc -> x.1.mrc. Based on A/pix and mass (in kd), the x-file will be scaled so an isosurface threshold of 1 will contain ~ the specified mass. 'ali to' is an iteration number. ie if 'ali to' is 4, then x.7.mrc would be aligned in 3D to x.4.mrc. x.3.mrc would not be aligned at all. Often this is set to a large value, like 99.
- 3dit=<iter>, 3dit2=<iter> No longer suggested. They apply a real-space iterative reconstruction technique to the 3D model to clean up artifacts caused by Fourier techniques. Typically 3dit=1 and 3dit2=2. Rarely necessary, and can produce other artifacts.
- amask=<r>,<threshold>,<iter> This option applies an automatically generated 'form fitting' soft (Gaussian) mask to the model after each iteration The mask generation is generally quite good. See proc3d option automask2 for details. This option can only be used in conjunction with xfiles=, since selection of the threshold requires proper volume normalization. This option can have a profound effect on proper convergence, but should be used with caution. Too tight a mask will produce artificially inflated resolution values. To use properly, insure that the mask does not cut through any significant density in the final model, and that the mean value is already ~0 at the mask position.

TYPICAL (not required):

- shrink=<n> Another option that can • produce dramatic speed improvements. In some cases, this option can actually produce an improvement in classification accuracy. This option scales the particles and references down by a factor of n before classification. Since data is often heavily oversampled, and classification is dominated by low resolution terms, this can be both safe, and actually improve classification by 'filtering' out high resolution noise. Generally shrink=2 is safe and effective especially for early refinement. In cases of extreme oversampling, larger values may be ok. This option should not be used for the final rounds of refinement at high resolution.
- euler2=<oversmp factor> This option should produce improvements in convergence and reconstruction quality, but has produced mixed results in the past. Not recently tested. It adds an additional step to the refinement process in which class-averages orientations are redetermined by projection-matching. The parameter allows you to decrease the angular step (ang=) used to generate projections. ie - 2 would produce projections with angular step of ang/2. It may be worth trying, but use it with caution on new projects. Untested with some new refine options.

EXPERIMENTAL (not required):

- usefilt This flag allows one to use arbitrarily filtered raw particles for classification purposes, but still use the unfiltered data when generating the actual reconstruction. To use this option, apply filter the data from start.hed into start.filt.hed. This option is now recommended for most reconstructions where ctfcw= is used. The particles should be filtered with proc2d wiener=.
- collapse=<alt max> This option is used to correct for a problem wherein side views of particles with Dn symmetry can iteratively 'drift' from alt=90 up to smaller altitudes. This problem should no longer exist if dfilt, phasecls or fscls are used.
- perturb This is a very new option, which is potentially useful, and at worst should be harmless. However, it has not been well characterized yet. Rather than generating Euler angles at evenly spaced positions, it adds some randomness to the positions. This should produce a more uniform distribution of data in 3D Fourier space and reduce Fourier artifacts.
- imask=<rad> This option was designed to improve classification for virus particles containing non-icosahedral DNA/RNA. It has not been tested in a long time and may or may not function.